

ECO PHYSICS CLD 88

Application examples



- Ambient monitoring in areas with excellent air quality
- Supervision of production processes in the chemical and high technology industries
- Permanent monitoring of clean room conditions in R & D labs
- Biomedical and pharmaceutical research
- Plant physiological research

The CLD 88 nitrogen oxide analyzer is unique in its speed and precision. It makes possible the continuous measurement of NO concentrations even in the range of parts per trillion!



Clean room laboratories require reliable and precise gas analysis.

When decimals are decisive.

The CLD 88 fulfills the requirements of many research laboratories specializing in detecting and monitoring smallest variations of NO concentrations in less than a second despite its total sample flow. The lagtime of less than a second makes it even more attractive.

NO_x is measured by an optional molybdenum converter. Unwanted effects of long sampling lines are minimized by the electronic and mechanical bypass system (option r).

User friendliness.

The development of an ECO PHYSICS analyzer always includes full user comfort. The user can adapt the operation according to his needs and applications by selection of predefined settings.

Warning messages or malfunctions are displayed coded and in full text. The analyzer guides the user step by step to return to normal operation.

The use of first-rate components guarantees virtually service-free operation. Maintenance simply means annual replacement of filters and membranes besides the consumables required by special sampling conditions.

Unique calibration by pressing a button!

The accuracy of chemiluminescence detection is strongly dependent on the calibration of the analyzer.

In order to assure reliability of its results the CLD 88 analyzer has optionally a calibration module (I) for the zero level and the NO reference gas. Calibration is quickly and automatically carried out by pressing a button on the keypad. This extremely useful feature eliminates the potential risk of erroneous calibrations.

Compact and modular construction.

The CLD 88 is the most compact unit of its class. Thanks to the totally modular layout and the rich variety of options this analyzer is designed for a multitude of applications.

- Compact design without any additional space required
- Optional molybdenum converter for NO_x detection
- Four freely selectable measurement ranges
- Operation and control via keypad or personal computer
- Optional bypass system to increase the sample flow
- Rapid system integration
- Optional calibration module for zero level and span gas calibration.



```
NO Range SNGL
39.32 100.0 Ppb
```

```
Meas.Mode Recorder Autocal MENU
Date/Time Comm. Timers Filter
```

```
Time for Calibration: 60 [s] MENU
NO-Cylinder Conc.: 100.0 [ppm]
```

```
Stand-by [V/MERS] ?
Zenair Pressure out of Range W-02
```


Clearly structured and full text displays inform the user about the instrument's status, any errors and measures to return to normal operation.

CLD 88

Specifications

| | | | |
|--|---|--|---|
| Measuring ranges | four freely selectable ranges from 5–5000 ppb | Analog output | 4–20 mA into 500 Ω max.; 0–1 V; 0–10 V |
| Min. detectable concentration | 0.05 ppb* | Dimensions | height: 133 mm (5¼") width: 450 mm (19") with moulding: 495 mm depth: 545 mm |
| Noise at zero point (1 σ) | 0.025 ppb* | Weight | 24 kg |
| Lagtime | <1 sec | Delivery includes | CLD 88 analyzer, power cable, analog signal cable, manual |
| Rise time (0–90%) | <1 sec | Standard | CLD 88 NO analyzer |
| Temperature range | 5–40 °C | Options | Y molybdenum converter I automatic calibration module for zero level and span reference gas r electro-mechanical pressure regulation |
| Humidity tolerance | 5–95% rel. h (non-condensing, ambient air and sample gas) | * depending on filter setting | |
| Sample flow rate | 0.3 l/min (1.2 l/min with option r) | ECO PHYSICS reserves the right to change these specifications without notice. | |
| Input pressure | ambient | | |
| Dry air use for O ₃ generator | internally generated (no external supply gas required) | | |
| Power required | 400 VA (incl. membrane pump and ozone scrubber) | | |
| Supply voltage | 100 - 230 V / 50 - 60 Hz | | |
| Interface | RS 232 | | |

Flow diagram

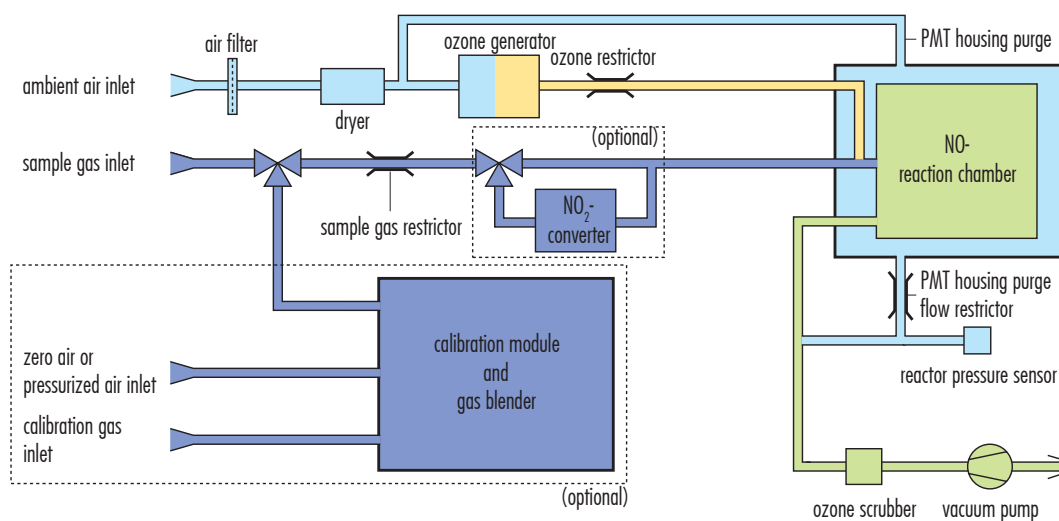

ECO PHYSICS

ECO PHYSICS INC. · 3915 Research Park Drive, Suite A-3 · ANN ARBOR, MI 48108-2200 · USA · Phone: (734) 998-1600

sales@ecophysics-us.com · www.ecophysics-us.com